Step 2: Installing Eyewash Unit

1. Remove Allen screw from Eyewash Yoke Assembly using the enclosed small Allen wrench (see Figure 3).

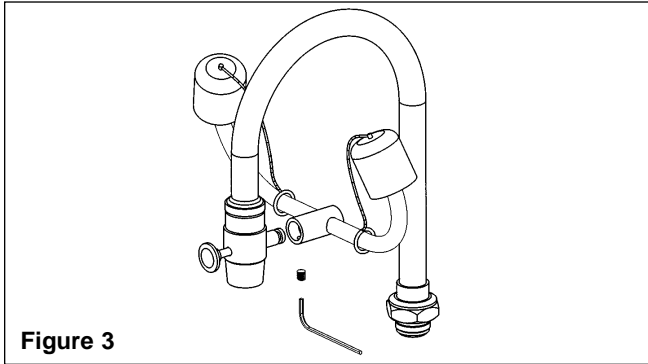
Figure 3

2. Attach Eyewash Yoke Assembly to connection on Eyewash adapter (see Figure 4). Be sure that Eyewash Spray Heads are facing toward the user's eyes.
3. Re-install Allen screw using small Allen wrench. Tighten Allen screw until eyewash yoke assembly is firmly attached to the eyewash adapter.

NOTE: When adjusting Eyewash Yoke assembly on adapter, always remember to loosen Allen screw first, make adjustment and then tighten.

4. Activate cold water supply and pull actuator pin to ensure eyewash is working properly. Yellow spray head dust caps will pop off with the pressure of the water. Make sure dust caps are on spray heads when eyewash is not in use.
5. Place Emergency Eyewash Sign and Inspection Tag in plain view by the eyewash.

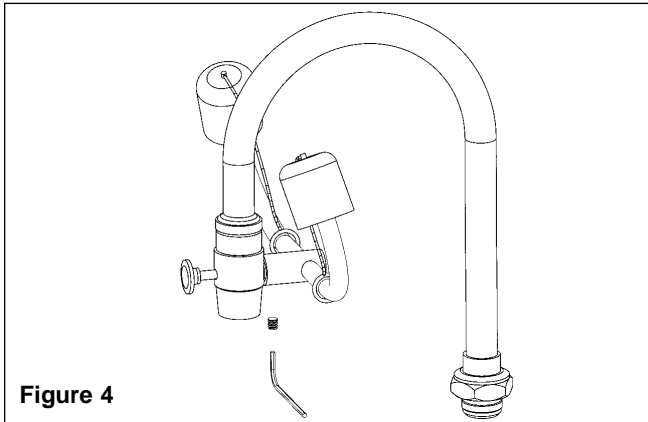
Figure 4

Step 3: Using Eyewash Unit

1. Turn cold or tempered water supply on.
2. Pull the actuator pin toward you.
3. Use the forefinger and thumb to hold your eyes open during rinse.
4. Continue to rinse eyes for a minimum of 15 minutes.
5. Seek medical attention immediately.

IMPORTANT


It is the responsibility of the employer to provide the following information to its employees:

- Circumstances that would render the use of an eyewash station.
- Location of each eyewash station in your facility.
- Proper use of the eyewash station, including the exclusive use of the cold water supply if tempered water cannot be provided (please refer to the Warning stated in the Preinstallation Information).

For further information on the proper installation and use of Emergency Eyewash and Shower equipment, please refer to ANSI Z358.1-2004.

**OBSOLETE
ECN 07-522**

Faucet-Mount Personal Eyewash Model S19-200

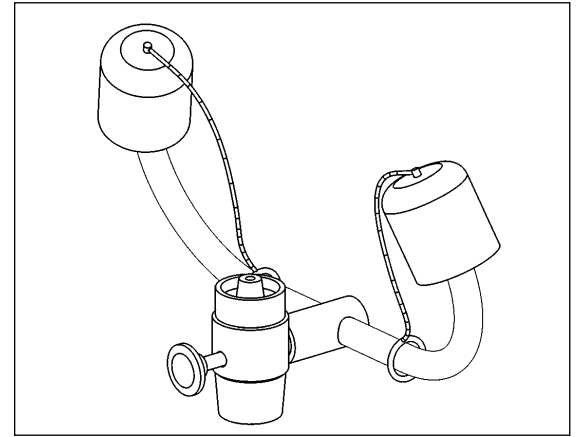Installation Instructions


CAUTION: The Faucet Mount Eyewash, S19-200 is considered a Personal Wash Unit, a type of supplemental equipment. Personal Wash Units do not meet the criteria of plumbed or self-contained eyewash equipment. Supplemental Equipment supports plumbed or self-contained units but shall not replace them.*

*ANSI Z358.1-2004

Bradley 

P.O. Box 309, Menomonee Falls, WI 53052-0309
TEL. 1-800-BRADLEY FAX 262-251-5817
<http://www.bradleycorp.com> Made in Taiwan

Model No. S19-200 Faucet-Mount Eyewash Parts List

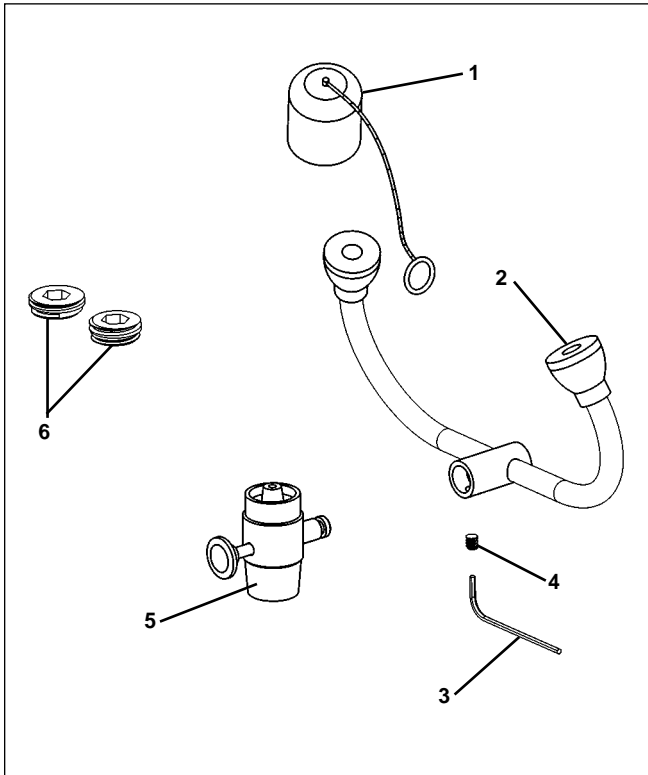Item	Part No.	Qty	Description
1	107-450	2	Dust Cover
2	S90-350	1	Yoke Assembly
3	130-140	1	Allen Wrench
4	160-131	1	Set Screw
5	S27-290	1	Diverter Assembly
6	S65-154	1	Threaded Adapter Kit
7*	S65-153	1	Gasket Kit
8*	204-486	1	Emergency Sign
9*	204-421	1	Inspection Tag

* Not shown

Pre-Installation Information

Check Contents

Separate all parts from the packaging and check each part with the parts breakdown on the preceding panel. Make sure all parts are accounted for before discarding any packaging material. If any parts are missing, do not attempt to install your Bradley Faucet-Mount Eyewash Unit until you obtain the missing parts.

Supplies Required:

- Wrench
- Standard or Gooseneck Faucet with 55/64"x 27 thread, 15/16"x 27 thread or 13/16"x 27 thread connections.

Note: Due to the many types of faucet designs available, the enclosed thread connections may not fit your faucet. It may be necessary to purchase the correct thread connections locally. Bradley has provided a 55/64"x27 thread connection on the Faucet Mount Eyewash adapter. Two additional thread connections (15/16"x27 and 13/16"x27 thread) have been provided to fit most common faucet designs. For additional adapter sizes, please contact Bradley.

- Cold or tempered water supply (minimum 30 PSIG)

WARNING: To prevent scalding of the eyes when eyewash is in use, it is recommended that hot water supply be disconnected OR that tempered water be provided. A comfortable eyewash water temperature range is 15° C - 35° C (60° F - 95° F). In circumstances where chemical reaction is accelerated by water temperature, a facilities safety/health advisor should be consulted regarding the optimum temperature for each application*.

*ANSI Z358.1-2004

Step 1: Installing adapter

1. Remove the existing aerator from the faucet (see Figure 1).

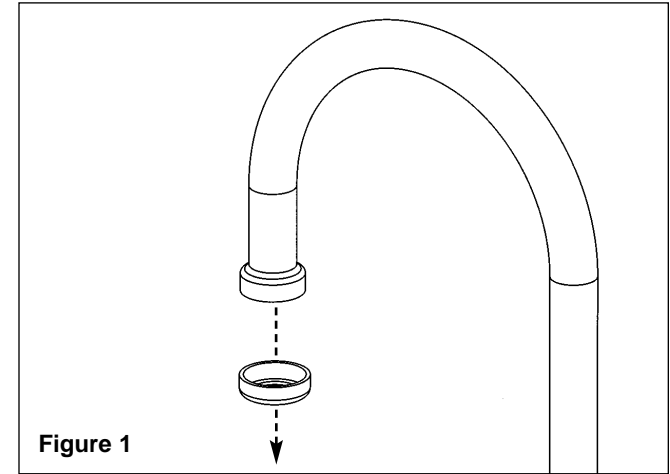
Figure 1

2. Screw the Faucet Mount Eyewash adapter onto the faucet and tighten with a wrench (see Figure 2). Actuator Pin (A) must be facing front, toward the user.

Note: To prevent scratches on the Eyewash adapter, it is recommended a grip pad or cloth be placed around the adapter while tightening with wrench.

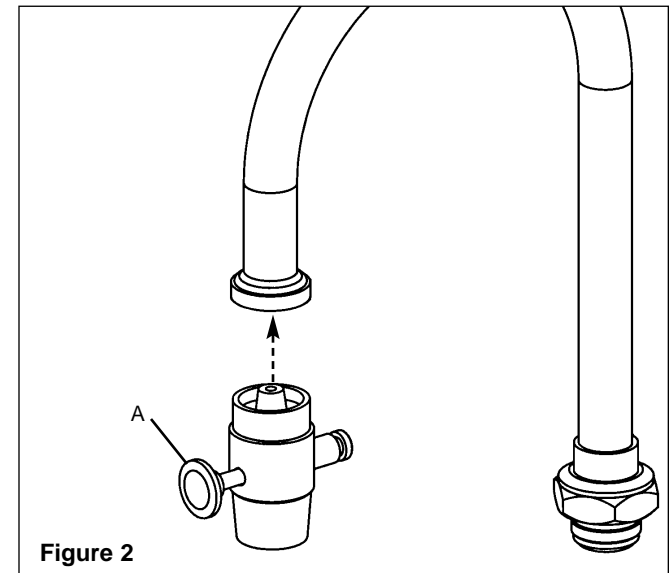
Figure 2

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