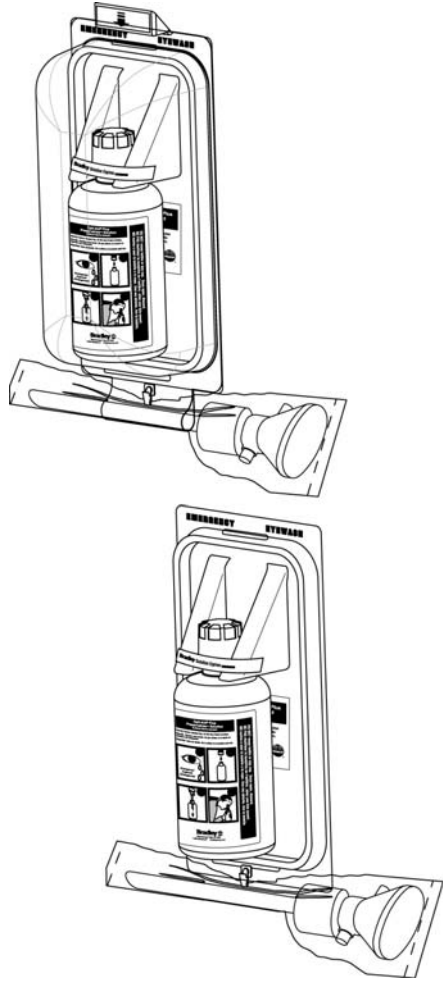
Installation

**Opti-Aid® Plus Personal Eyewash Station
with PreservaFree™ Solution**

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Table of Contents

Pre-Installation Information	2
Directions for Use	3
Installation	4-5

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**S19-945 (32 oz. Opti-Aid® Plus Refill) (Remplir)
(Recambio)**

⚠ WARNING ⚠

Infection, loss of vision or other serious injury may result from misuse. **DO NOT** use expired, cloudy, colored or contaminated solution. **DO NOT** use if safety seal is broken. **DO NOT** save unused solution. **DO NOT** refill bottle. Instruct employees of bottle location and how to properly use eyewash bottle in case of an emergency.

Personal eyewash equipment does not meet the criteria of plumbed or self-contained eyewash equipment. Personal eyewash equipment supports but shall not replace plumbed or self-contained eye-wash equipment. This product is intended to provide first aid relief prior to reaching a plumbed or self-contained eyewash station. Please refer to ANSI Standards or local area safety codes for specific eye and eye/face wash equipment requirements and their intended use.

For questions regarding the operation or installation of this product, visit www.bradleycorp.com or call 1-800-BRADLEY. Product warranties may also be found under "Product Information" on our web site at www.bradleycorp.com.

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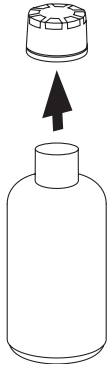
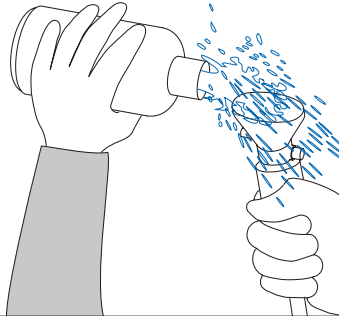
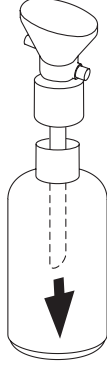
1

2

FOR UNITS WITHOUT DUST COVER: Secure the holder to the wall with two fasteners provided by installer. Proceed to Step 3.

FOR UNITS WITH DUST COVER: Secure the holder to the wall with two fasteners provided by installer. **NOTE: DO NOT tighten the bottom fastener at this time!** Proceed to Step 2.

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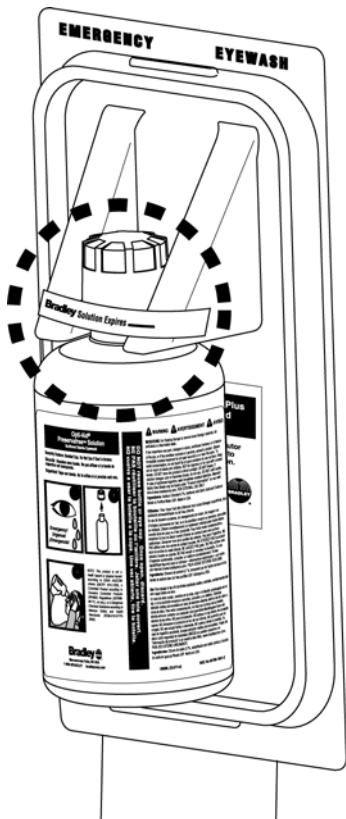
FOR UNITS WITH DUST COVER: Slide the dust cover hook behind the holder and down over the fastener. Tighten the bottom fastener, securing the dust cover to the wall.

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
3

4

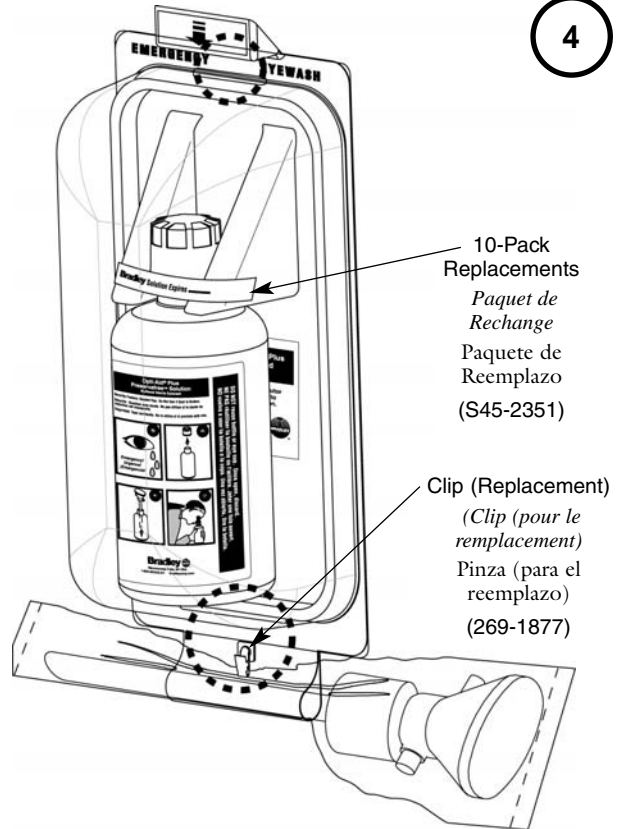FOR ALL UNITS: Position the bottle in the holder with “Opti-Aid Plus” instructions facing outward. Write the month and year your bottled eyewash solution expires onto the “Solution Expires” label. The expiration date is located on the neck of the bottle. Attach the “Solution Expires” label to the holder.

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FOR UNITS WITHOUT DUST COVER: Secure the clip to the holder, then secure the bag to holder with the clip.

FOR UNITS WITH DUST COVER: Secure the clip to the dust cover, then secure the bag to holder with the clip. Position the dust cover and press the upper snap to secure. While holding the dust cover in place, press the lower snap.

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